
## KIMTECH PURE\* A5 Sterile Cleanroom Apparel with CLEAN-DON\* Technology



- **Step 1:** (Pre-Entry) Don Hair Net and Shoe Covers after removing all jewelry and cosmetics.
- Step 2: (Gowning) Wash hands and gown first pair of sterile gloves. Sanitize gloves after gowning each article if required.
- Step 3: Apply mask and hood assuring a snug fit.



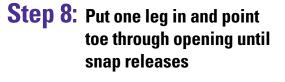

**Step 9:** Do the same with the other leg

**Step 12:** Slip thumbs through thumb loops



**Final Step:** Cross legs and zip up coverall

