

KIMTECH PURE* A5 Sterile Boot Cover with Grasp Ties Donning Technique



Step 1

Select your boot cover size; inspect package for vacuum seal; tear along notches toward the product to open; flex package to separate boot covers



Step 2

Pull out one boot cover by the cuff



Step 3

Grasp ties with one hand and make sure they don't touch the floor



Step 4

Open up boot cover; point toe towards opening and pull up over calf while holding the ties



Step 5

Hold up foot or rest on bench. Grasp ties at front of boot cover and wrap over cusp of shoe/boot; then wrap ties behind ankle and tie in front.



Step 6

Place donned boot on the clean side



Step 7

Pull the remaining boot cover out of package by the cuff and repeat Steps 3, 4 and 5. Proceed with final gowning steps



Exceptional
Workplaces*

